


# Infinity XE 50X1 Series

User Manual

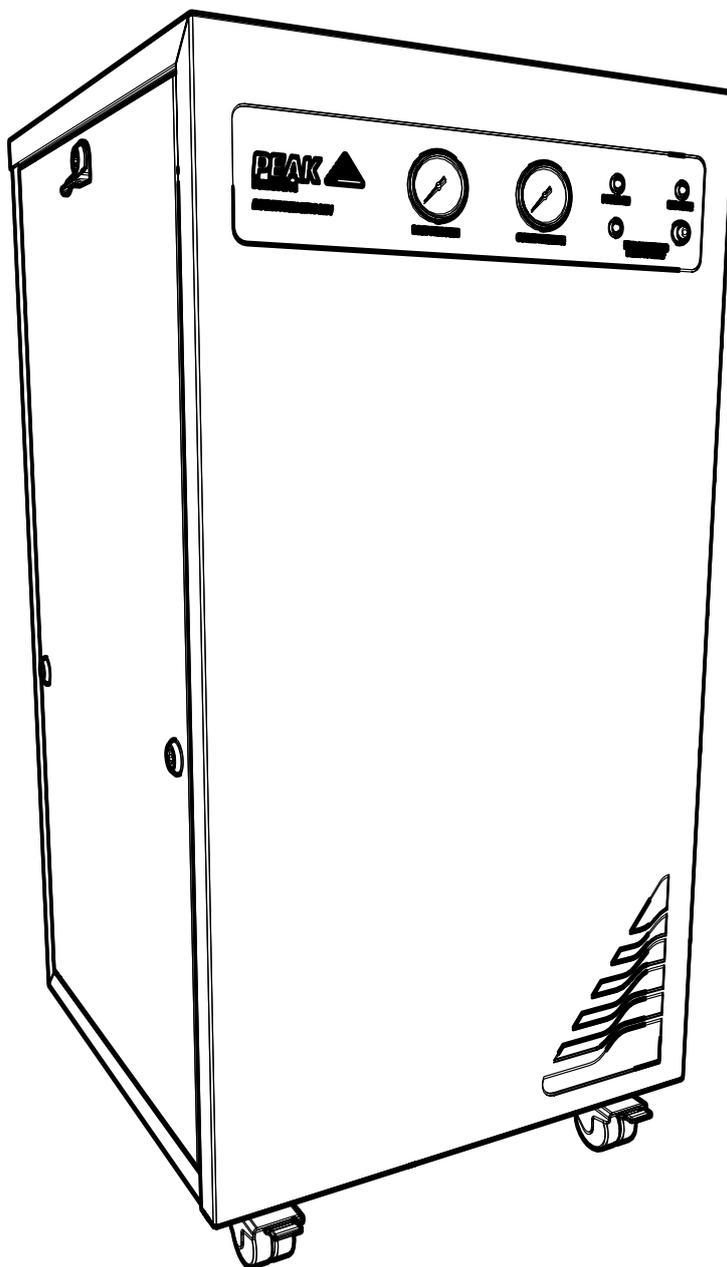

# Contents

Change History	3
How to use this Manual	3
Introduction	4
Warranties and Liabilities	5
Warranty & Liability Coverage	5
Safety Notices	7
Symbols	7
Safety Notice to Users	7
Declaration of Conformity	8
Environmental Declaration	9
Technical Specification	10
Unpacking	12
Fittings Kit Contents	13
Installation	14
Generator Environment	14
Generator Overview	15
General Dimensions	15
Rear Connections	16
Inlet Air Quality	16
Electrical Connection	17
Electrical Panel Access	17
Normal Operation	18
On Demand Gas	18
O <sub>2</sub> Alarm	18
Eco Mode	18
Commissioning	19
Oxygen Analyser	19
Tubing Lengths	19
Service Requirements	20
Service Schedule	20
Peak Protected	21
Cleaning	22
Safely Isolated Condition	22
Safe State After Repair	22
Troubleshooting	23

## Change History

Rev	Comment	Name	Date
2	Infinity XE 5031 & 5041 Changes	Liam Couttie	29/11/2017

## How to use this Manual

This manual is intended for end users and has been written as a reference document where you can skip to the relevant information.

Users can refer to the contents page to find the relevant information.

Please review each of the following sections carefully.

Thank you for selecting Peak Scientific to meet your gas generation needs, and should you require any further assistance or support please do not hesitate to contact Peak Scientific or the Peak Partner from which you purchased your generator.

# Introduction

The Infinity XE 50 Series delivers performance and value beyond expectations, offering a pure and reliable source of gas. These perpetual flow nitrogen generators are suitable for a variety of applications in your laboratory.

These models provide a source of nitrogen gas with other features including:

- Durable - Few moving parts with minimum maintenance requirements
- Single Source Solution - One generator can cater for the requirement of multiple applications & multiple laboratories on site
- Quiet in operation - Completely silent whilst operating
- Mobile - Flexibility to position the generator where required
- Simple installation - Generator designed as a plug and play system
- Economical - More cost effective than any other gas supply method
- Convenient - Gas on demand, no hazards, no need to worry about running out of gas
- Oxygen Analyser - Built-in purity monitoring system with safe guard technology
- Eco-mode - Energy efficient eco-mode ensures low running costs

The Infinity XE 50 Series generators are based on proven technology, and selectively remove oxygen, moisture and other gases to leave clean, dry, phthalate free nitrogen.

To ensure these Generator models meet our high expectations with regards to reliability and performance, we have tested them extensively at our manufacturing plant and with end users around the world to ensure reliability and longevity of the system.

# Warranties and Liabilities

## Warranty & Liability Coverage

1. Peak warrants that, subject to the provisions in this statement, purchased Peak generators, whether purchased directly from Peak or indirectly via an approved, certified and trained distributor or partner (referred to hereafter as a "Peak Partner") will comply in all material respects with any specifications referred to in your customer order confirmation and, subject to installation and operational guidelines being followed as described in applicable product manuals, shall be free from any defects in quality of materials or workmanship for a period of one year from the date of installation, provided this takes place within 3 months of factory dispatch.
2. Where the purchased generator is from the Precision Hydrogen series, Peak further warrants that, subject to installation and operational guidelines being followed as described in applicable product manuals, the hydrogen cell shall be free from any defects in quality of materials or workmanship for a total period of three years (inclusive of warranty period specified in clause 1) from date of installation, provided this takes place within 3 months of factory dispatch.
3. Where the purchased generator is from the i-Flow 6000 series, Peak further warrants that, subject to installation and operational guidelines being followed as described in applicable product manuals, the generator shall be free from any defects in quality of materials or workmanship for a total period of two years (inclusive of warranty period specified in clause 1) from the date of installation, provided this takes place within 3 months of factory dispatch and the following provisions have also been met:
  - a. you must purchase a service plan, ensuring the generator is serviced by Peak or a Peak Partner on or before the end of the first 12 months of your ownership, and serviced at least once during each subsequent 12 month period thereafter;
  - b. the generator (and any associated equipment) must have been commissioned by Peak or a Peak Partner;
  - c. the feed air or inlet air supply to the generator must comply with ISO 8573-1:2010 Class 1.2.1 at all times;
  - d. your air compressor, dryer, filtration and oil removal systems must be deemed suitable for use by Peak or a Peak Partner, and must be changed and serviced regularly, in line with the equipment manufacturer's recommended guidelines; and
  - e. any generator failure or fault that is deemed to have been caused by the failure of any upstream equipment, component, part or system (such as air compressor, air treatment or filtration) will be excluded from the warranty described herein.
4. Peak also warrants that any replacement parts whether purchased (directly from Peak, or via a Peak Partner) or supplied as part of any remedial action undertaken in line with the provisions of clauses 12 and 13, shall be free from any defects in quality of materials or workmanship for a period of 180 days from the date of factory dispatch, provided its installation is performed by Peak or a Peak Partner.
5. This warranty does not exclude Peak's liability in respect of any claim for death or personal injury to any person, in so far as such can be attributed to negligence or breach of duty of care directly resulting from failure of Peak to comply with the provisions in clauses 1, 2, 3 & 4.

## Exclusions & Limitations

6. This warranty does not cover:
  - a. damage, deterioration or malfunction resulting from an alteration or modification to a generator which has not been carried out by Peak or a Peak Partner;
  - b. damage, deterioration or malfunction resulting from what Peak reasonably believes to be abuse, or misuse of a generator by you or any third party;
  - c. liability for accident or neglect (other than pursuant to clause 5);
  - d. maintenance or repairs which have not been carried out by Peak or a Peak Partner;
  - e. operation of a generator or exposure of a generator to environmental conditions that fall out-with operational guidelines as specified in the applicable product user manual; and
  - f. lightning, power surges or any other acts of God or nature.
7. This warranty is non-transferrable. Only the original owner of the generator may benefit from the terms within this statement.

8. Peak shall not be liable in respect of any claim made for costs, damages, losses or expenses (whether consequential, direct, indirect or otherwise) or in any respect howsoever arising including, but not limited to, liability from accident or negligence (other than pursuant to clause 5) that may be suffered by you or any third party.
9. No person or entity is authorised to change the terms and conditions outlined in this warranty statement in any respect, or to create any additional obligations or liabilities for any party involved.
10. This warranty statement supersedes any and all prior warranty agreements between the parties and constitutes the complete, final and exclusive understanding of the parties with respect to the subject matter. All prior negotiations, representations, or promises, whether oral or written, of either party shall be deemed to have been merged herein.
11. If any part of this warranty statement is invalidated, for whatever reason, such part will be deleted and the rest shall remain unaffected, continuing to be in full force and effect.

### **Delivery of Warranty Service**

12. Subject to clause 13, and:
  - a. Peak being notified by you, within the duration of the applicable warranty period, of any defect that you think is subject to any warranty valid under clauses 1, 2, 3 or 4; and
  - b. Peak being permitted to inspect the generators, parts and their installation (along with any relevant packaging)  
Peak shall at its option repair or replace defective generators or parts (including, if necessary, any moving parts and irrespective of runtime). No additional charges will apply, for parts or delivery and, where applicable, labour or travel. Peak will endeavour to deliver this service within 3 working days of your notification.
13. Where, in Peak's reasonable opinion, a defect is subject to an exclusion described in clause 6, Peak reserves the right to charge for parts or delivery and, where applicable, you may also be charged by Peak for call out, labour or travel in respect of any repair or replacement which you authorize Peak to carry out.

# Safety Notices

Peak Scientific Instruments cannot anticipate every possible circumstance which may represent a potential hazard. The warnings detailed within this manual refer to the most likely potential hazards, but by definition cannot be all inclusive. If the user employs an operating procedure, item of equipment or a method of working which is not specifically recommended by Peak Scientific, the user must ensure that the equipment will not be damaged or become hazardous to persons or property.

## Symbols

This manual uses the following symbols to highlight specific areas important to the safe and proper use of the generator.

 WARNING	A WARNING notice denotes a hazard. It calls attention to an operating procedure, process or similar, which if not correctly performed or adhered to, could cause personal injury or in the worst case death. Do not proceed beyond a WARNING notice until the indicated conditions are fully understood or met.
 CAUTION	A CAUTION notice denotes a hazard. It calls attention to an operating procedure, process or similar, which if not correctly performed or adhered to, could cause damage to the generator or the application. Do not proceed beyond a CAUTION notice until the indicated conditions are fully understood or met.
	Caution, risk of electric shock. Ensure power to the generator has been removed before proceeding.

## Safety Notice to Users



These instructions must be read thoroughly and understood before installation and operation of your Peak Generator. Use of the generator in a manner not specified by Peak Scientific MAY impair the SAFETY provided by the equipment.



When handling, operating or carrying out any maintenance, personnel must employ safe engineering practices and observe all relevant local health and safety requirements and regulations. The attention of UK users is drawn to the Health and Safety at Work Act 1974, and the Institute of Electrical Engineers regulations.



If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment maybe impaired.

# Declaration of Conformity

We Peak Scientific Instruments Ltd.  
Of Fountain Crescent, Inchinnan, Renfrewshire, PA4 9RE

Declare that:

Equipment: Nitrogen Gas Generator  
Models: Infinity XE 50 Series

To which this declaration relates, is in conformity with the applicable EC directives, harmonized standards, and other normative requirements.

- Low Voltage Directive 2014/35/EU  
EN 61010-1: 2010 Electrical Equipment for measurement, control and laboratory use.
- Electromagnetic Compatibility Directive 2014/30/EU  
EN 61326-1: 2013 Electrical Equipment for measurement, control and laboratory use.

All evaluation, testing and certification issued by:

## York EMC Services

Unit 1, Grangemouth Technology Park  
Earls Road  
Grangemouth  
UK  
FK3 8UZ

Signed:  Name: Chris Pugh

Date: 22/06/2017 Position: Chief Technology Officer



# Environmental Declaration

We Peak Scientific Instruments Ltd.  
Of Fountain Crescent, Inchinnan, Renfrewshire, PA4 9RE

Declare that:

Equipment: Nitrogen Gas Generator

Models: Infinity XE 50 Series

Is fully compliant with the following Directives

- 2012/19/EU WEEE (Waste of Electrical and Electronic Equipment)
- 2011/65/EU RoHS 2 (Restriction of Hazardous Substance)

Peak Scientific Instruments Ltd fully complies with its obligations towards the European WEEE (Waste of Electrical and Electronic Equipment) Directive 2012/19/EU. These obligations are being met within the B2B compliance group.

Peak Scientific Instruments Ltd has developed all reasonable 'due diligence' controls to ensure that our products comply with the principles and requirements of the European recast RoHS (Restriction of Hazardous Substances) Directive 2011/65/EU. Similar directives in the United States and China, for example, have also been captured within this program.

Where a specific certificate of compliance is required, this can be requested, on a product serial number basis, directly from Peak Scientific Instruments Ltd, by contacting us through our website on [www.peakscientific.com](http://www.peakscientific.com)

Signed:  Name: Chris Pugh

Date: 22/06/2017 Position: Chief Technology Officer



# Technical Specification

## Infinity XE 5011 & 5021

### Environment

	5011	5021
Minimum Operating Ambient Temperature	5°C (41°F)	
Maximum Operating Ambient Temperature	30°C (86°F)	
Maximum Relative Humidity	80%	

\* Maximum safety ambient temperature 30°C

\*\* When taken out of storage the generator should be allowed to acclimatize at room temperature for a minimum of 3 hours before operation.

### Inlet Conditions

Minimum Air Inlet Pressure	4.1 bar (60 psi)
Maximum Air Inlet Pressure	9.9 bar (145 psi)
Minimum Air Inlet Flow	Dependant on required flow and purity
Minimum Air Quality	ISO 8573-1:2010 Class [1.4.1]

### Generator Outlets

Maximum Pressure Drop (Inlet-Outlet)	0.7bar (10psi)	
Maximum Outlet Flow (95% Nitrogen)	130 L/min	260 L/min
Outlet Nitrogen Purity	95 - 99.5%	
Start-Up Time	30 minutes	
Particles	<0.01µm	
Pressure Gauges	2	

### Electrical Requirements

Voltage	100-230V	
Frequency	60 Hz / 50 Hz	
Current	0.4 A / 0.174 A	
Input Connection	C20 Plug	
Power Cord (Supplied)	C19 socket to local connection (13A minimum)	
Pollution Degree	2	
Insulation Category	I	

### General

Dimensions cm (inches) WxDxH	50 x 50 x 99 (19.6 x 19.6 x 39.0)	
Generator Weight Kg (lbs)	50.4 kg (111 lb)	54.6 kg (120 lb)
Shipping Weight Kg (lbs)	67 kg (147 lb)	72 kg (159 lb)

# Technical Specification

## Infinity XE 5031 & 5041

### Environment

	5031	5041
Minimum Operating Ambient Temperature	5°C (41°F)	
Maximum Operating Ambient Temperature	30°C (86°F)	
Maximum Relative Humidity	80%	

\* Maximum safety ambient temperature 30°C

\*\* When taken out of storage the generator should be allowed to acclimatize at room temperature for a minimum of 3 hours before operation.

### Inlet Conditions

Minimum Air Inlet Pressure	4.1 bar (60 psi)
Maximum Air Inlet Pressure	9.9 bar (145 psi)
Minimum Air Inlet Flow	Dependant on required flow and purity
Minimum Air Quality	ISO 8573-1:2010 Class [1.4.1]

### Generator Outlets

Maximum Pressure Drop (Inlet-Outlet)	1.4bar (20psi)	
Maximum Outlet Flow (95% Nitrogen)	390 L/min	520 L/min
Outlet Nitrogen Purity	95 - 99.5%	
Start-Up Time	30 minutes	
Particles	<0.01µm	
Pressure Gauges	2	

### Electrical Requirements

Voltage	100-230V	
Frequency	60 Hz / 50 Hz	
Current	0.4 A / 0.174 A	
Input Connection	C20 Plug	
Power Cord (Supplied)	C19 socket to local connection (13A minimum)	
Pollution Degree	2	
Insulation Category	I	


### General

Dimensions cm (inches) WxDxH	50 x 50 x 99 (19.6 x 19.6 x 39.0)	
Generator Weight Kg (lbs)	58.8 kg (130 lb)	77 kg (170 lb)
Shipping Weight Kg (lbs)	63 kg (139 lb)	82 kg (181 lb)

# Unpacking

Although Peak Scientific takes every precaution with safe transit and packaging, it is advisable to fully inspect the unit for any sign of transit damage.

Check 'SHOCKWATCH' and 'TIP-N-TELL' labels for signs of rough handling prior to unpacking

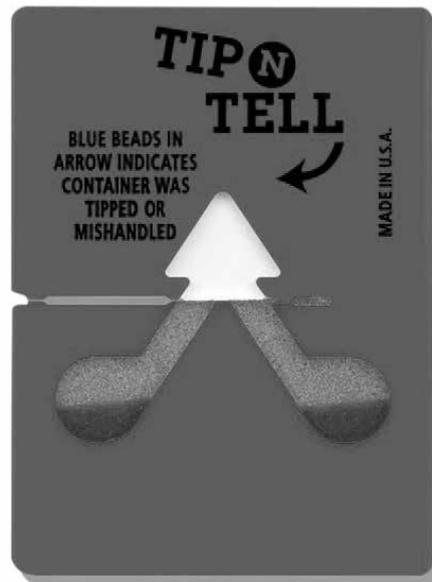

**Any damage should be reported immediately to the carrier and Peak Scientific or the Peak Partner from where the unit was purchased.**

Follow the unpacking instructions posted on the side of the crate. It will require two people to remove the unit from the shipping crate and to manoeuvre the generator to the desired location.

Please save the product packaging for storage or future shipment of the generator.

Note: Included with the generator is a "Fittings Kit" containing mains power leads for UK, EU & US and also all the required fittings and warranty registration card. Be careful not to discard these with the packaging.

## Fittings Kit Contents

Supplied in the Fittings Kit are all the fittings required to connect the generator to the application. The contents of the Fittings Kit are as follows:

1. 6mm PE Tubing	x 3m
2. 12mm PE Tubing	x 6m
3. 6mm Push-Fit Fitting	x 1
4. 12mm Push-Fit Fitting	x 2*
5. 16mm Push-Fit Fitting	x 1**
6. Tee fitting	x 1
7. Shut-off Valve	x 1
8. UK Mains Power Cable	x 1
9. EU Mains Power Cable	x 1
10. JP Mains Power Cable	x 1
11. US Mains Power Cable 230v	x 1
12. US Mains Power Cable 110v	x 1
13. 4mm Hex Key	x 1
14. 5mm Hex Key	x 1
15. 8mm Hex Key	x 1

All of the generators output ports are located on the output panel at the rear of the unit.

\* x1 with 5031 & 5041

\*\* 5031 and 5041 only

# Installation

## Generator Environment

The generator is designed for indoor use only. It should be installed adjacent to the application(s) it is supplying. If this is not convenient then the unit can be sited elsewhere, however, consideration should be made of the lengths of pipe runs as pressure drops can result from extended runs of pipe.


Performance of the generator (like all sophisticated equipment) is affected by ambient conditions. Note should also be taken to the proximity of Air Conditioning outlets. These can sometimes give rise to “pockets” of air with high relative humidity. Operation of the unit within such a pocket could adversely affect its performance. Consideration should also be given to the air flow around the unit. It is recommended that an air gap of 75mm (3”) should be maintained down both sides and at the rear of the unit. Please refer to the drawing below for the general dimensions of the unit.

Minimum Operating Ambient Temperature:        5 °C (41 °F)

Maximum Operating Ambient Temperature:        30 °C (86 °F)


# Generator Overview

## General Dimensions

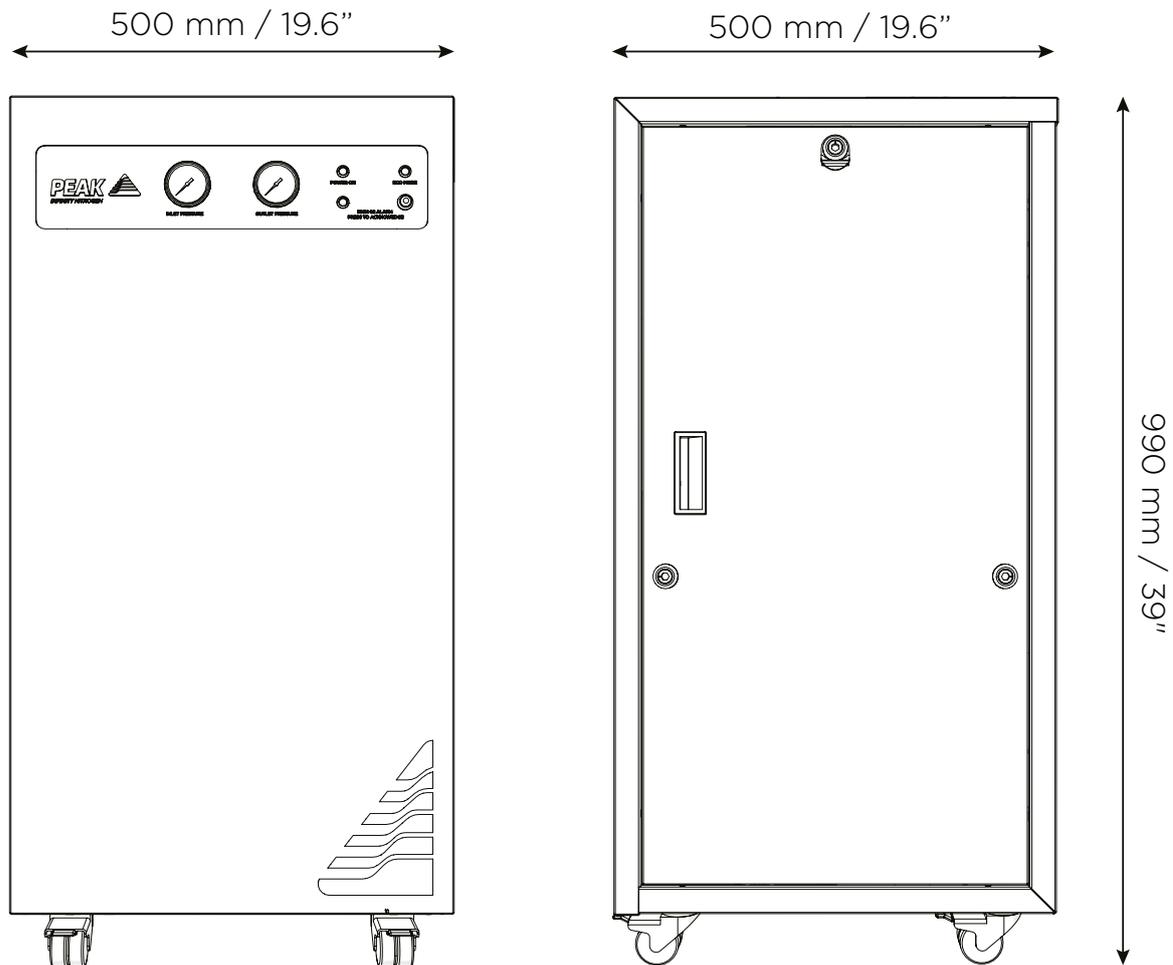

The generator must always be placed on a flat, level surface. Failure to do so will affect the performance of the generator.

## Rear Connections

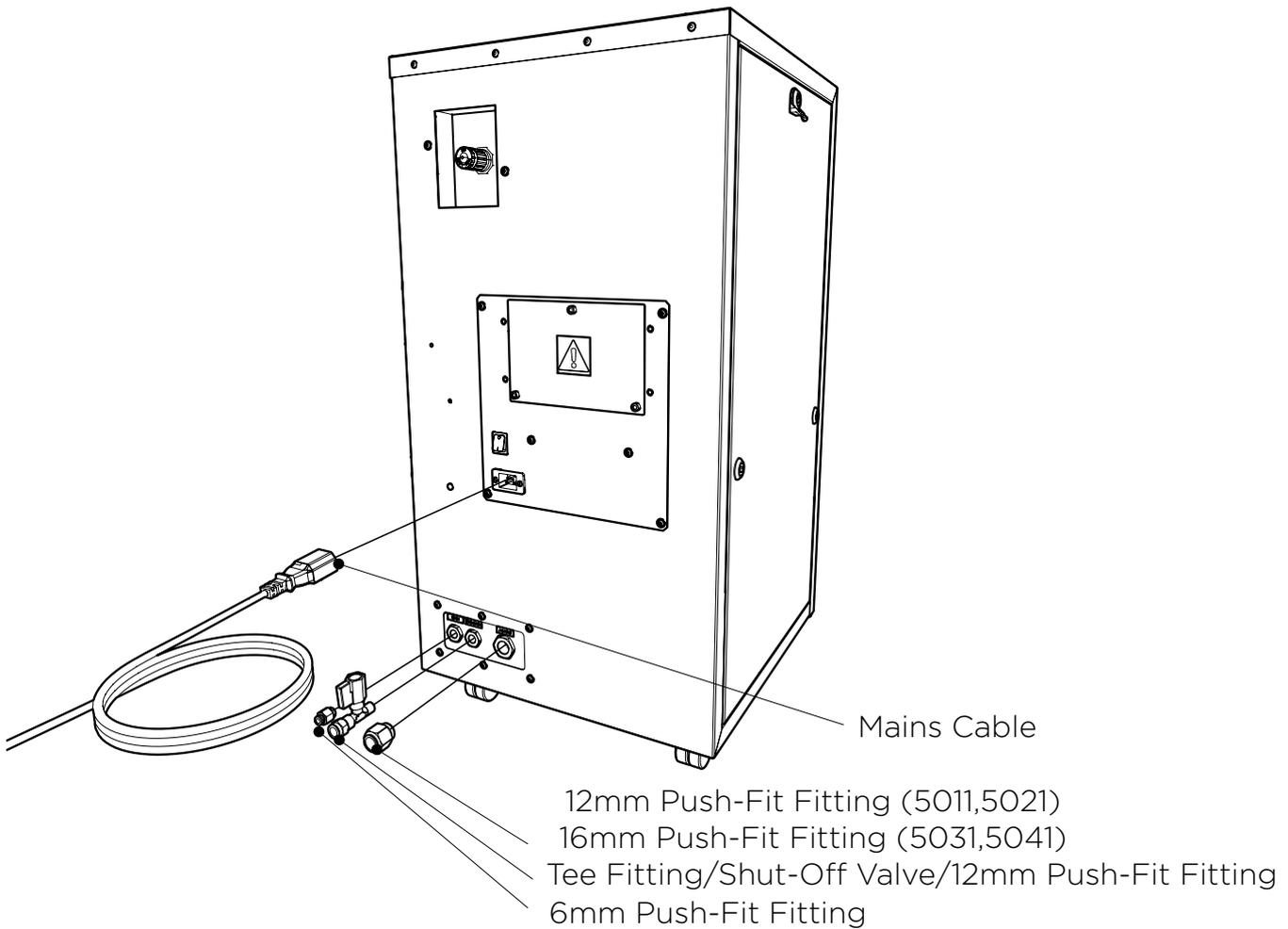

## Inlet Air Quality

The Nitrogen Generator should be connected to an ISO 8573-1:2010 (1:4:1) source of compressed air. A minimum pressure of 60 psig is required for efficient operation of the Generator. Any doubts as to the suitability of your compressed air supply should be referred to the factory for advice.

### **A MAXIMUM OIL VAPOUR CONTENT OF 0.01 mg per m<sup>3</sup> is PERMISSIBLE**

Air Inlet and Nitrogen Outlet connections are at the rear of the generator in the middle & right side of the panel. The Air inlet connection is the right port, 12mm Push-Fit Fitting (5011,5021) 16mm Push-Fit Fitting (5031,5041), Outlet port is 12mm Push-Fit Fitting

# Electrical Connection

Connect the generator to an appropriate 85-264 volt single-phase supply, refer to the generator serial plate for input specification and ensure your supply matches the requirements.

If the appropriate power cord is not supplied; a new plug, rated to at least 12 amps, can be fitted by a qualified electrician.



This unit is classified as SAFETY CLASS 1. THIS UNIT MUST BE EARTHED. Before connecting the unit to the mains supply, please check the information on the serial plate. The mains supply must be of the stated AC voltage and frequency.

EARTH/GROUND (E):-                      Green & Yellow      or      Green

LIVE (L):-                                      Brown                      or      Black

Neutral (N):-                                  Blue                                  or      White

# Electrical Panel Access



Changing settings within this enclosure may adversely affect the operation of the generator. This enclosure should not be accessed unless by or under direct instruction from a trained individual.

## Normal Operation

The Infinity XE 50 Series Gas Generators are designed specifically to minimize operator involvement. Given that the systems are installed as described in earlier sections and is serviced in accordance with the specified maintenance recommendations (see Service Requirements), then it should simply be a matter of turning the Generator on when it is required.

The Generator will automatically produce the factory set flow and pressure as detailed in the Technical Specifications.

## On Demand Gas

The generator will produce nitrogen on demand. If the application is operating and requires a gas flow, the system it is connected to will supply this to suit the requirements of the application. If the application requirement for gas stops, the system will also stop. If the demand from the application starts again, the system will detect the demand for gas and will automatically start again to suit the demand.

## O<sub>2</sub> Alarm

The O<sub>2</sub> analyser live monitors output gas purity against a user-specified limit. If the oxygen concentration exceeds this limit, the generator's oxygen warning light will be illuminated. If oxygen concentration exceeds the specified limit for 30 minutes, the generator will close down and depressurise to protect purity-sensitive applications. For generators supplying pressure-sensitive applications, this auto-shutdown can be overridden to ensure applications always see the required pressure.

## Eco Mode

When demand from the instrument is below 1.2L/min the Infinity XE generator will stop demanding air from the external compressor. If demand from the application starts again, the system will detect the demand for gas and will automatically start up again.

## Commissioning

With the generator installed as previously described, connect the application to the Nitrogen output port and open the air supply to the unit. On reaching operational pressure the generator will produce Nitrogen at the customer-specified purity. The design of the generator is such that it will deliver up to the rated output flow of Nitrogen, determined by demand of the consuming equipment. Rated flow, purity and pressure are customer-specific and dependent on inlet conditions being met as detailed on the individual generator

The design of the generator is that it will deliver variable purity dependant on the inlet pressure and outlet flow rate. Outlet pressure may be adjusted by the pressure regulator on the rear of the unit, allowing for up to 15 psi pressure drop across the system. Should the demand for the nitrogen stop, the generator will automatically go into "eco-mode" (refer to page 18), shutting off the air inlet. Once the demand restarts the generator automaticcally resumes producing Nitrogen. The generator is protected from over pressurising.

## Oxygen Analyser

The generator has the option to activate an Oxygen Analyser that can be set to activate a warning if the oxygen level is higher than the specified level. If the light is illuminated this means that the Oxygen level is higher than the defined threshold. If the Oxygen level is consistently high for more than 30 minutes and the reset button has not been pressed, the generator will automatically shut-down. The reset button will return the unit to normal operation. If the oxygen level remains high, the alarm and shut down will repeat.

## Tubing Lengths



The diameter of the tubing which will be connected to the gas outlet is important and is determined by the length of tubing required. Failure to follow these recommendations could lead to pressure between generator and application.

> 10 - 40 meters: Use 12/10 (12mm O/D, 10mm I/D) Tubing.

> 40 metres: Please contact Peak Scientific with the relevant distance and we will calculate the flow resistance and the tubing size required.

The imperial equivalentents are: 12/10 = 1/2" O/D, 3/8" I/D.

**Air input tubing for 5031 & 5041 is not supplied. 16mm OD, 12mm ID PTFE Tubing is recommended.**

# Service Requirements


## Service Schedule

Purchase Interval	Component	Visit
12 Months	Infinity XE 50 Series Annual Maintenance Kit	<a href="http://www.peakscientific.com/ordering">www.peakscientific.com/ordering</a>

# Peak Protected

With Peak Scientific you invest in not only a product but peace of mind. With a network of certified Peak engineers stationed throughout the globe, Peak's rapid response team are never far away and our commitment is to keep your generator running day in, day out, protecting your laboratory workflow.

**[Peak Protected]** can provide...

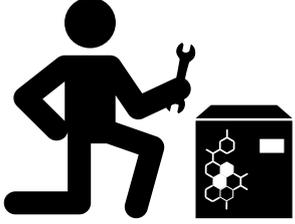


**Installation**  
A dedicated Peak engineer will visit your lab to install and setup your generator




**Complete plan**  
Swift response by a Peak Service Engineer within 72 hours & planned preventative maintenance



**Premium Protected**  
Guaranteed rapid on-site response within 24 hours & planned preventative maintenance

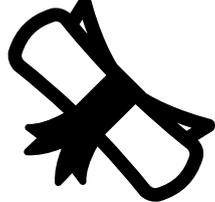

**IQ/OQ**  
Certified assurance for applications requiring documented qualification

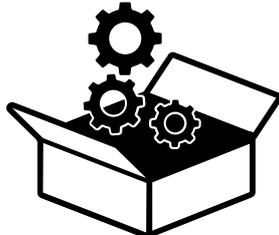

**Spares**  
Genuine Peak parts with express delivery, ensuring optimal performance and lifetime



**Technical Support Hotline**  
Around the clock support by phone or online with our global technical helpdesk

To find out more about protecting your investment visit: [www.peakscientific.com/protected](http://www.peakscientific.com/protected)

## Cleaning

Clean the outside of the generator only using warm soapy water and a clean damp cloth. Ensure all excess fluid is thoroughly removed from the cloth prior to use.



Cleaning should only be undertaken with the power switched off and the power cord removed from the rear of the generator.



Under no circumstances should any solvents or abrasive cleaning solutions be used as these can contain fumes that could be harmful to the generator.



Care should be taken with Leak Detections Liquids.

## Safely Isolated Condition

The unit is in a safely isolated condition when it is disconnected from its application and fully de-pressurised. Directions for isolating the Generator are shown below.



Failure to place the Generator in a safely isolated condition when instructed to do so may lead to personal injury or injury to others and even death.

1. Switch generator off.
2. Disconnect from air supply.
3. Ensure the output pressure gauge reads zero. (If gauge does not fall to zero, open the manual ball valve at the outlet port, to allow trapped gas to escape.)
4. Disconnect from the application.

## Safe State After Repair

To guarantee the generator is in a safe state after a repair, please ensure the following conditions are met;

- All earth leads have been reconnected.  
Earth leads should be connected to the earth tabs located on the door panels, as shown below.
- The electricity cable has been correctly refitted.  
The mains cable should be plugged into the IEC power inlet located on the rear panel of the generator. **This unit is classified as SAFETY CLASS 1. THIS UNIT MUST BE EARTHED.**

# Troubleshooting

Problem	Possible Solution
<p>The Generator will not switch on and the power switch does not illuminate.</p>	<ul style="list-style-type: none"> <li>• Ensure power cord is plugged into the Generator and that the power socket is turned on.</li> <li>• Check the fuse in the power cord plug.</li> <li>• Contact your service provider.</li> </ul> 
<p>The Generator will not switch on but the power switch is illuminated.</p>	<ul style="list-style-type: none"> <li>• Disconnect power cord from the rear of the Generator. Open the left hand panel are check that both circuit breakers are turned on (switch in the up position). Reconnect power cord.</li> <li>• Contact your service provider.</li> </ul> 
<p>The output pressure has dropped below its original set point.</p>	<ul style="list-style-type: none"> <li>• Check the pressure regulator at the back of the unit has not been adjusted.</li> <li>• Check to see if the inlet pressure gauge has dropped below the set air supply pressure.</li> <li>• If so, check your air supply does not have a pressure drop over the length of tubing and has no leaks.</li> <li>• Contact your service provider.</li> </ul>
<p>The generator will not switch on but the power indicator and O2 alarm are illuminated.</p>	<ul style="list-style-type: none"> <li>• Disconnect applications from the generator to reduce demand to within rated capacity.</li> <li>• Press O2 alarm acknowledgement button.</li> <li>• Contact your service provider.</li> </ul>

# Go Online or Complete and Return

We know that registering any of your recently purchased products is not the first thing on your mind- but it is very important to both of us. Not all warranties are alike and Peak Scientific stand out against other gas suppliers as we offer a comprehensive, quick response, on-site warranty. This means that in the very unlikely case that your gas generator develops a fault we have rapid support teams on-hand around the world who are able to come to your lab and get you back up and running in no time.

Register for your **comprehensive 12 month on-site warranty** with ease online at [www.peakscientific.com/protected](http://www.peakscientific.com/protected).

Alternatively, you can send the completed form to Peak Scientific by post or email at [warranty@peakscientific.com](mailto:warranty@peakscientific.com).

**Go Online or Complete and Return**

You can register for your **FREE 12 month Warranty** with ease online at [www.peakscientific.com/protected](http://www.peakscientific.com/protected).

Alternatively, you can send the completed form to Peak Scientific by post or email at [warranty@peakscientific.com](mailto:warranty@peakscientific.com).

Product Warranty Registration			
Contact name			
Email address			
Company			
Address			
City/town			
Postcode			
Country			
Telephone			
Generator serial #			
Model type			
Installation date			
Do you still use an alternative gas solution i.e. cylinders or bulk liquid?	Yes	No	
What gas requirements do you have in your lab?	Hydrogen	Nitrogen	Zero Air


**Extend your cover with**

Peak Scientific offer comprehensive gas generator after sales support packages. Peak [Protected] aftercare support can guarantee an on-site response within 72 hours\*, genuine parts from our ISO9001 approved factory and a 95% first-time fix rate. See our enclosed Peak [Protected] leaflet for further information.

**Important!**

You have 1 month to register your Peak Scientific product from the date of installation. Once registered the warranty will be honoured for a period of 12 months. If you wish to defer the installation of your generator, you must notify Peak Scientific immediately by emailing [warranty@peakscientific.com](mailto:warranty@peakscientific.com). For generators that remain unregistered after 1 month from the shipment date, the warranty will be considered active from the date of factory dispatch.

\* Complete Plan only

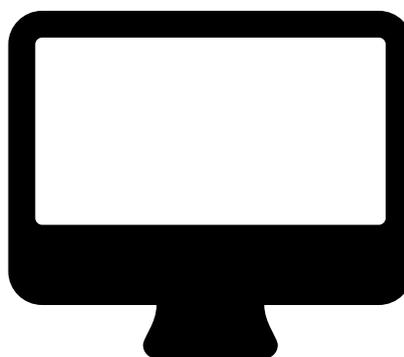

## Important!

You have **1 month to register** your Peak Scientific product from the date of installation. Once registered the warranty will be honoured for a period of 12 months. If you wish to defer the installation of your generator, you must notify Peak Scientific immediately by emailing [warranty@peakscientific.com](mailto:warranty@peakscientific.com). For generators that remain unregistered after 1 month from the shipment date, the warranty will be considered active from the date of factory dispatch.







# [**PEAK Protected**]<sup>TM</sup>

Peak Scientific has highly trained, fully certified Field Service Engineers located in over 20 countries across every continent around the world. This allows us to provide an industry-leading rapid response service to our customers. With **[Peak Protected]**, your laboratory's productivity becomes our top priority.

To discuss Peak Protected generator cover and payment options speak to your local Peak Representative or for further information contact: [protected@peakscientific.com](mailto:protected@peakscientific.com)

## **Peak Scientific**

Fountain Crescent  
Inchinnan Business Park  
Inchinnan  
PA4 9RE  
Scotland, UK  
**Tel:** +44 141 812 8100

For further information on any of our generator products please contact [marketing@peakscientific.com](mailto:marketing@peakscientific.com)

