

# Genius 3040 User Manual



## Contents

Change History	4
How to use this Manual	4
Introduction	5
Warranties and Liabilities	6
Safety Notices	7
<i>Symbols</i>	7
Safety Notice to Users	7
EU Declaration of Conformity	8
UK Declaration of Conformity	9
WEEE Compliance Statement	10
Technical Specification	11
Environment	11
Generator Outlets	11
Electrical Requirements	11
General	11
Unpacking	12
Installation	13
Generator Environment	13
Removal of Transit Brackets	14
Unit Controls	15
Rear Connections	15
Fittings Kit	16
Drain Connection	16
Electrical Connection	17
Start-up Sequence	19
Connecting to the Mass Spectrometer	20
Tubing Lengths	20
IMPORTANT DOCUMENTS	21
Normal Operation	22
On Demand Gas	22
Generator Cycling	22

Unusual Operation	22
Service Requirements	23
Service Schedule	23
Service Indication	23
Stage 1	23
Stage 2	24
Service Indication Reset	24
Service Plans	24
Cleaning	24
High Duty Indication	25
High Duty Indication Reset	25
Indication of Fault	25
Trouble Shooting	26

## Change History

Rev.	Comment	Name	Date
1	Change to Drain Port Location	Liam Couttie	14/06/2013
2	Capacitor Change	Liam Couttie	18/03/2014
3A	Content Update	Liam Couttie	30/03/2015
4	Fittings Kit Update	Liam Couttie	05/02/2016
5	Declarations Update	Liam Couttie	16/11/2021

## How to use this Manual

This manual is intended for end users and has been written so that it can either be read as a step by step guide to installation and usage or as a reference document where you can skip to the relevant information.

Users of a hard copy version can refer to the contents page to find the relevant information. Users of the soft copy version can use the hyperlinks from the contents page as well as the hyperlinks between sections.

Please review each of the following sections carefully.

Thank you for selecting Peak Scientific to meet your Gas Generation needs, and should you require any further assistance or support please do not hesitate to contact Peak Scientific or Peak Partner from which you purchased your Generator.

## *Introduction*

The Genius 3040 has been developed to cater for the requirements of the SCIEX range of LCMS instruments.

This model provides two independent sources of Nitrogen gas as well as two independent supplies of clean, dry air; other features include:

Other features include:

- Quiet in operation – noise level of 54 db(A)
- Anti- vibration – maximum reduction of vibration
- Service indication – allowing you to plan your maintenance and keep your LCMS uptime at a maximum
- Improved drainage – reduction of moisture carry over and thus increased reliability
- Re- heat technology – improves membrane performance and reliability
- Robust control system – improves safety and reliability of units

With the Genius 3040 based on proven technology, it selectively removes oxygen, moisture and other gases to leave clean, dry, phthalate free Nitrogen and Dry Air. Internal air compressors make this unit independent from in- house air supplies and fitted castors allow the user to easily position the unit in the lab. The unit has further been approved by SCIEX to supply curtain, source and exhaust gas to all of their LCMS applications.

To ensure this Generator model meets our high expectations with regards to reliability and performance, we have tested this new model extensively at our manufacturing plant and with end users around the world to ensure reliability and longevity of the system.




## *Warranties and Liabilities*

1. The Company warrants that it has title to the Goods.
2. Subject to the provisions of this clause the Company warrants that the Goods shall comply in all material respects with any specification referred to in the Order Confirmation (as the same may be amended) and shall, subject thereto, be free from defects in material and workmanship for the lesser of a period of twelve months from the date of delivery or thirteen months from the date of dispatch from the factory.
3. Save as provided in this clause and except where the Goods are sold to a person dealing as a consumer (within the meaning of the Unfair Contract Terms Act 1977) all warranties, conditions or other terms implied by statute or common law are hereby expressly excluded save to the extent they may not be lawfully excluded. When the Goods are sold to a consumer within the meaning of the Unfair Contract Terms Act 1977 their statutory rights are not affected by the provisions of this clause.
4. In the event of the Customer making a claim in respect of any defect in terms of clause 2 hereof the Customer must.
  1. Reasonably satisfy the Company that the Goods have been properly installed, commissioned, stored, serviced and used and without prejudice to the generality of the foregoing that any defect is not the direct or indirect result of lack of repair and/or servicing, incorrect repair and/or servicing, use of wrong materials and/or incorrect spare parts
  2. Allow the company to inspect the Goods and/or any installation and any relevant packaging as and when reasonably required by the Company.
5. Subject to the Company being notified of any defect as is referred to in sub-clause 2 hereof within a reasonable time of it becoming apparent and subject always to the terms of sub-clause 4 hereof, the Company shall, in its option, replace or repair the defective Goods or refund a proportionate part of the Price. The Company shall have no further liability to the Customer (save as mentioned in sub-clause 6 hereof).
6. The Company shall be liable to indemnify the Customer in respect of any claim for death or personal injury to any person in so far as such is attributable to the negligence or breach of duty of the Company or any failure by the Company to comply with the provisions of sub-clause 2 hereof.
7. Save as provided in sub-clause 2 hereof the Company shall not be liable in respect of any claim by the Customer for costs, damages, loss or expenses (whether direct, indirect, consequential or otherwise) or indemnity in any respect howsoever arising including, but not by way of limitation, liability arising in negligence (other than pursuant to clause 6 above) that may be suffered by the Customer or any third party.

## Safety Notices

### Symbols

This manual uses the following symbols to highlight specific areas important to the safe and proper use of the Generator

	A <b>WARNING</b> notice denotes a hazard. It calls attention to an operating procedure, process or similar, which if not correctly performed or adhered to, could cause personal injury or in the worst case death. Do not proceed beyond a <b>WARNING</b> notice until the indicated conditions are fully understood or met.
	A <b>CAUTION</b> notice denotes a hazard. It calls attention to an operating procedure, process or similar, which if not correctly performed or adhered to, could cause damage to the Generator or the Application. Do not proceed beyond a <b>CAUTION</b> notice until the indicated conditions are fully understood or met.
	Caution, risk of electric shock. Ensure power to the Generator has been removed before proceeding.

### Safety Notice to Users



*These instructions must be read thoroughly and understood before installation and operation of your Peak Genius 3040 Generator. Use of the Generator in a manner not specified by Peak Scientific MAY impair the SAFETY provided by the equipment.*



*When handling, operating or carrying out any maintenance, personnel must employ safe engineering practices and observe all relevant local health and safety requirements and regulations. The attention of UK users is drawn to the Health and Safety at Work Act 1974, and the Institute of Electrical Engineers regulations.*

## EU Declaration of Conformity

We Peak Scientific Instruments Ltd.  
Of Fountain Crescent, Inchinnan, Renfrewshire, PA4 9RE

Hereby declare that, this declaration of conformity is issued under the sole responsibility of the manufacturer.

Equipment Type: Nitrogen & Dry Air Generator  
Model Designator: Genius 3040

To which this declaration relates, is in conformity with the following applicable EU Directives, harmonized standards, and other normative requirements.

- **Low Voltage Directive 2014/35/EU**  
EN 61010-1: 2010 Safety Requirements for Electrical Equipment for Measurement, Control and Laboratory Use.
- **Electromagnetic Compatibility Directive 2014/30/EU**  
EN 61326-1: 2013 Electrical Equipment for Measurement, Control and Laboratory Use - EMC Requirements. (Class A)
- **Restriction on the use of certain hazardous substances in electronic equipment (RoHS) Directive 2011/65/EU as amended by EU 2015/863.**

Signed for and on behalf of Peak Scientific by

Signed:



Name: Fraser Dunn

Position: Design Engineering Manager  
Peak Scientific Instruments Ltd,  
Inchinnan, Renfrew, Scotland, PA4 9RE, UK.

Date: 10th August 2021





## UK Declaration of Conformity

We Peak Scientific Instruments Ltd.  
Of Fountain Crescent, Inchinnan, Renfrewshire, PA4 9RE


Hereby declare that, this declaration of conformity is issued under the sole responsibility of the manufacturer.

Equipment Type: Nitrogen & Dry Air Generator  
Model Designator: Genius 3040

To which this declaration relates, is in conformity with the following applicable UK Statutory Instruments, Standards and other normative requirements.

- **The Electrical Equipment (Safety) Regulations 2016 (SI 2016 / 1101) as amended.**  
BS61010-1:2010 Safety Requirements for Electrical Equipment for Measurement Control and Laboratory Use.
- **The Electromagnetic Compatibility Regulations 2016 (SI 2016 / 1091) as amended.**  
BS61326-1:2013 Electrical Equipment for Measurement , Control and Laboratory Use - EMC Requirements.
- **The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012 (SI 2012 / 3032) as amended.**

Signed for and on behalf of Peak Scientific by

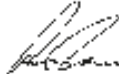Signed: 

Name: Fraser Dunn

Position: Design Engineering Manager  
Peak Scientific Instruments Ltd,  
Inchinnan, Renfrew, Scotland, PA4 9RE, UK.

Date: 10th August 2021



## WEEE Compliance Statement

The Waste Electrical and Electronic Equipment (WEEE) Regulations SI 2013 No 3113 and or the Waste Electrical and Electronic Equipment (WEEE) Directive 2012/19/EU apply to all electrical and electronic equipment placed on the market in the UK and EU covered by the scope of regulations which can be found in the Government Guidance Notes (PDF) produced by the Department for Business Innovation and skills for the UK and here for Europe.

All PEAK products that are subject to the WEEE directive are compliant with the WEEE marking requirement. Such products are marked with the "crossed-out wheelie bin" symbol (shown below) in accordance with European standard EN50419. All old electrical equipment can be recycled. Please do not dispose of any electrical equipment (including those marked with this symbol) in general rubbish bins. Please contact your dealer or distributor for clarity.



## Technical Specification

### Environment

Minimum operating ambient temperature	5°C (41°F)
Maximum operating ambient temperature*	35°C (95°F)
Maximum relative humidity	80% Non-Condensing
Maximum operating altitude	2000 meters
Minimum storage temperature**	-20°C (-4°F)
Maximum storage temperature**	60°C (140°F)

\* Maximum safety ambient temperature 35°C

\*\* When taken out of storage the Generator should be allowed to acclimatize at room temperature for a minimum of 3 hours before operation.

### Generator Outlets

Curtain maximum flow	12 L/min @ 5.5 bar (0.42 cfm @ 80 psi)
Source maximum flow	24 L/min @ 7.6 bar (0.84 cfm @ 110 psi)
Exhaust maximum flow	8 L/min @ 4.2 bar (0.28 cfm @ 60 psi)
Nitrogen maximum flow	32 L/min @ 6.8 bar (1.13cfm @ 100 psi)
Particles	<0.01µm
Phthalates	NONE
Suspended liquids	NONE
Gas outlets	4 x 1/4" BSPP
Drain outlet	1 x 1/4" BSPP
Pressure gauges	4

### Electrical Requirements

Voltage	230 VAC ± 10%
Frequency	50/60 Hz
Current	12 Amps
Input connection	C20 Plug
Power cord (Supplied)	C19 socket to local connection (13A minimum)
Pollution degree	2
Installation category	II

### General

Dimensions in cm (inches) W x D x H	60 x 85 x 133 (23.6 x 33.5 x 52.4)
Weight	199kg (439 lb)
Shipping weight	284 kg (626 lb)
Noise level	54 dBA @1m



*It should be noted that the gas pressures and flows are factory set. The pressures shown on the Generator front panel are in excess of the maximum inlet pressure of the Mass Spectrometer. This is to allow for known pressure drops. These settings have been approved by SCIEX.*

## Unpacking

Although Peak Scientific takes every precaution with safe transit and packaging, it is advisable to fully inspect the unit for any sign of transit damage.

Check 'SHOCKWATCH' label for signs of rough handling prior to un-packing –



*Any damage should be reported immediately to the carrier and Peak Scientific or the Peak Partner from where the unit was purchased.*

Follow the unpacking instructions posted on the side of the crate. It will require two people to remove the unit from the shipping crate and to manoeuvre the Generator onto the floor.

Please save the product packaging for storage or future shipment of the Generator.

Note: Included with the Generator is a "Fittings Kit" containing mains power leads for UK, EU and US also all the required fittings. Be careful not to discard these with the packaging.

## Installation

### Generator Environment




The Generator is designed for indoor use only. It should be installed adjacent to the Mass Spectrometer(s) it is supplying. If this is not convenient then the unit can be sited elsewhere, however, consideration should be made of the lengths of pipe runs as pressure drops can result from extended runs of pipe. Please see the [Tubing lengths](#) section for further details.



Performance of the Generator (like all sophisticated equipment) is affected by ambient conditions. Note should also be taken to the proximity of Air Conditioning outlets. These can sometimes give rise to “pockets” of air with high relative humidity. Operation of the unit within such a pocket could adversely affect its performance. Consideration should also be given to the air flow around the unit. It is recommended that an air gap of 75mm (3”) should be maintained down both sides, at the rear and across the top of the unit. Please refer to the drawing below for the general dimensions of the unit.

Maximum Ambient Conditions: 35°C (dry bulb) 80%RH (Max) Non-Condensing

### General Dimensions

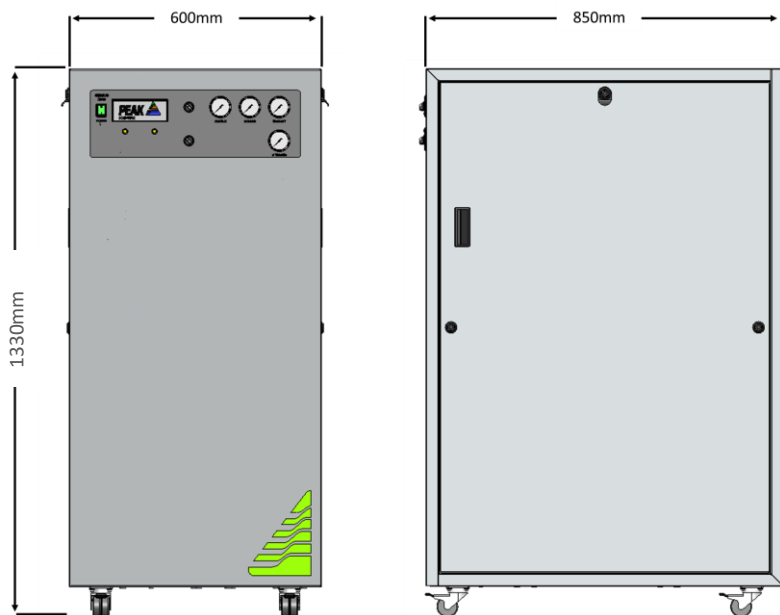

The Generator must always be placed on a level surface. Failure to do so will affect the performance of the Generator.

## Removal of Transit Brackets



*The transit brackets must be removed prior to switching the unit on. Failure to do so will result in damage to the equipment. This will void the warranty on the Generator and will result in a chargeable repair.*

1. Using the 8mm hex key from the Fittings Kit, remove the side covers from the cabinet ensuring that earth cables are disconnected.
2. Remove the four screws with a #3 Pozidrive Screwdriver. See Figure 1.
3. Slide the transit bracket out from under the compressor by pulling it towards you. See Figure 1.
4. Repeat this for the compressor on the other side of the Generator.
5. Retain the transit brackets as these must be refitted if the Generator is to be transported again.

Note. Do not re-fit door panels in preparation for Voltage Check

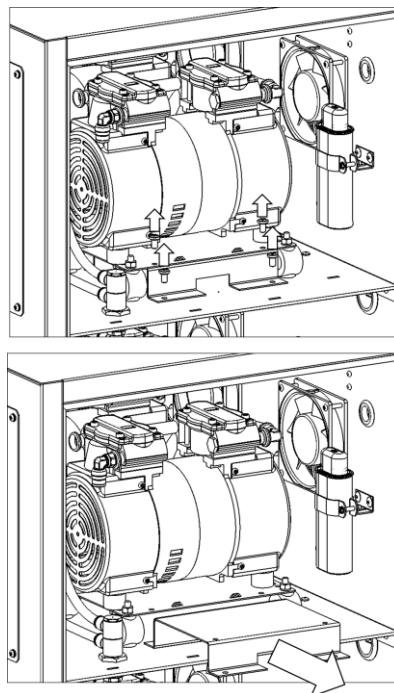


Figure 1: Transit Bracket Removal

## Unit Controls

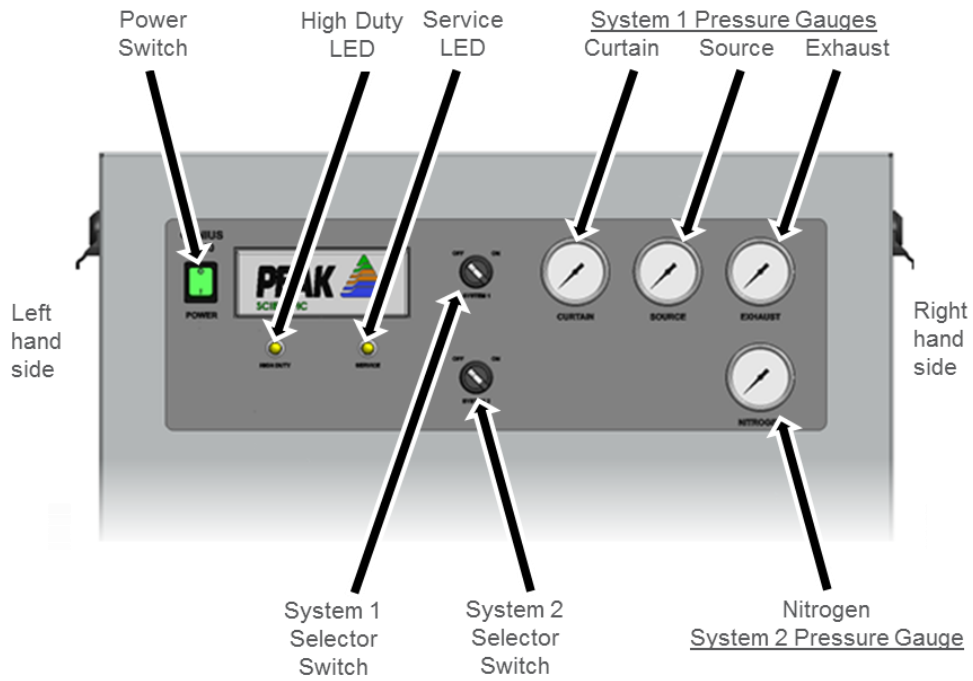


Figure 2: Unit controls

## Rear Connections

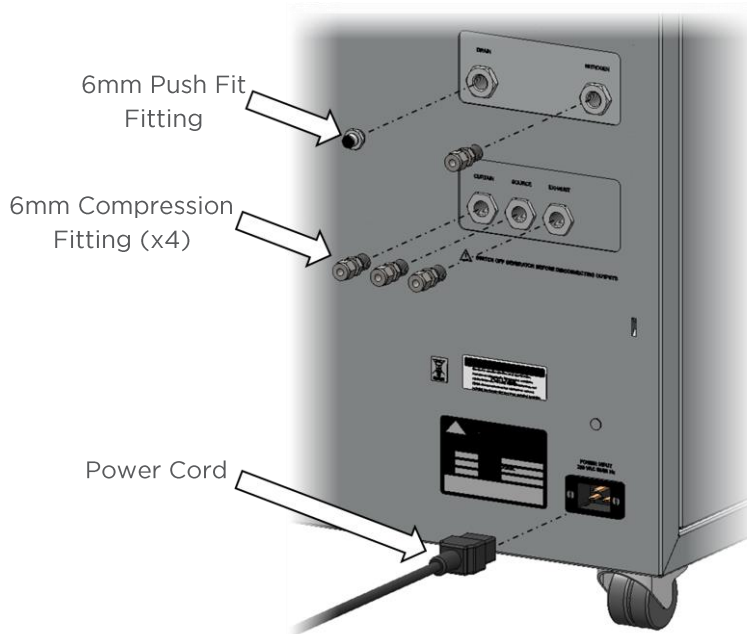


Figure 3: Rear Connections

## Fittings Kit

Supplied in the Fittings Kit are all the fittings required to connect the Genius 3040 Generator to the Mass Spectrometer(s). The contents of the Fittings Kit are as follows:

- |                                     |       |
|-------------------------------------|-------|
| 1. 8mm Hex Key                      | × 1   |
| 2. 6mm Push Fit Tube Fitting        | × 1   |
| 3. ¼" Compression Tube Fitting      | × 4   |
| 4. 6mm High Temp Tubing             | × 3m  |
| 5. ¼" Teflon Tubing                 | × 12m |
| 6. C19 Mains Cable (UK)             | × 1   |
| 7. C19 Mains Cable (Euro)           | × 1   |
| 8. C19 Mains Cable (US)             | × 1   |
| 9. Installation Guide – Genius 3040 | × 1   |

All of the Generator output ports are located on the output panel at the rear of the Generator.

## Drain Connection

Fit the 6mm push fit fitting (item 2 from fittings kit) to the drain port located on the output panel (see Figure 3). Tighten using a 16mm or 5/8" spanner. Use the 6mm tubing (item 4 from fittings kit) to connect this to a suitable drain connection or container. It should be noted that the Generator can expel a considerable amount of water from this (dependant on ambient humidity).



*If a container is used it should be emptied at regular intervals. The container must NOT have an air tight seal as water and air are expelled at pressure.*



## Electrical Connection

Connect the Generator to a 230 volt single-phase supply using the power cord provided. If the appropriate power cord is not supplied; a new plug, rated to at least 13 amps, can be fitted by a qualified electrician.



*This unit is classified as SAFETY CLASS 1. THIS UNIT MUST BE EARTHED. Before connecting the unit to the mains supply, please check the information on the serial plate. The mains supply must be of the stated AC voltage and frequency.*

EARTH/GROUND (E):-	Green & Yellow	or	Green
LIVE (L):-	Brown	or	Black
Neutral (N):-	Blue	or	White

Our electrical requirements are 230VAC nominal +/- 10%. This means that the Generator can accommodate transients between 207VAC and 253VAC. However, running continuously at voltages less than 220V is not recommended and extended periods at these extremes can have a detrimental effect on the operation and life of the Generator.

To ensure that the correct voltage is being supplied to the Generator, the Genius 3040 comes equipped with an inbuilt voltmeter. This is located on the underside of the compressor compartment, Figure 4 below, and will measure and display the mains voltage that is being supplied to the Generator.

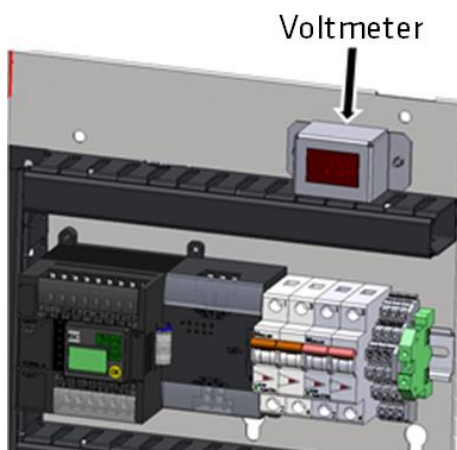


Figure 4: Voltmeter Location

The voltmeter should be checked prior to the initial purge of the system. To do this, with the doors remaining off and with the Generator still connected to the 230 volt single-phase supply, the unit should be powered ON using the power switch (identified by "I").



*Do not touch anything inside the Generator whilst the side panels are removed and the mains power is connected to the unit*

When the Generator is switched on, the input voltage will be displayed on the voltmeter. If this reading is 219V or less, then we would highly recommend fitting a transformer. This can be ordered directly from Peak Scientific with the order number being as below.

<i>Product Description</i>	<i>Part Number</i>
Dual Tap Transformer 200V - 230V	06-3210

*Table 1: Additional Transformer*

On completing the voltage check, the Generator should be powered OFF using the switch on the front panel, the power cord removed from the rear of the generator and the door panels refitted.



CAUTION

*Ensure the Earth wires are reconnected to the side covers when refitting.*

## Start-up Sequence

Once re-connected to the mains supply, the Generator can now be switched on. Flick the power switch on the front panel of the Generator to the on position (identified by “I”).



*Before the Generator is connected to the Mass Spectrometer the Generator should be operated in isolation (i.e. not connected to the Mass Spectrometer) for thirty minutes. This is to ensure any impurities present are purged from the system. Failure to do this may harm the Mass Spectrometer.*

Pressure will start to build in the internal storage tanks. This can be monitored by watching the relevant output pressure gauges on the front panel. These will climb to the factory set pressures as noted in the specifications.

Once these pressures are reached, the compressors will continue to run until the internal tank upper pressure limits (9.1 and 8.6 bar for Nitrogen and 8.9 bar for Dry Air) are reached and the compressor has run for a period of at least two and a half minutes. It should be noted that the Nitrogen and Dry Air compressors work completely separately from one another, thus their individual compressors may not have the same run periods.

The compressors will then rest until the internal tank lower pressure limits (7.9 bar for Nitrogen and 8.1 bar for Dry Air) are reached. Once this limit is reached the compressors will switch back on again. This compressor cycling is normal and will continue throughout the operation of the Generator.

When the system has been operated for a period of 30 minutes, all the internal pipe-work and storage tanks, will have been purged with clean Dry Air and Nitrogen.

The system is now purged and the tubes can be connected at the rear of the unit to the relevant Mass Spectrometer.



*Should a system be disconnected from its Mass Spectrometer, the 30 minute purge run must be performed before re-connecting the system to its specified Mass Spectrometer*

## Connecting to the Mass Spectrometer

Once the initial purge run of 30 minutes has completed, attach the 1/4" compression fittings supplied in the fitting kit to the Curtain, Source and Exhaust outlets on the output panel of the operating system. Using the Teflon tubing supplied connect each outlet of the Generator to the corresponding inlet on the Mass Spectrometer.

If you require more tubing than is supplied please refer to the [Tubing Lengths](#) section.



*It should be noted that the gas pressures and flows are factory set. The pressures shown on the Generator front panel are in excess of the maximum inlet pressure of the Mass Spectrometer. This is to allow for known pressure drops. These settings have been approved by SCIEX.*



*Once the tubing is connected to the Mass Spectrometer, please ensure that they are thoroughly checked for being leak-tight. Even the slightest leak in the gas supply between the Generator and the Mass Spectrometer can lead to a reduction in efficiency, or possible alarm of the Generator.*

## Tubing Lengths



*The diameter of the tubing which will be connected to the gas outlet is important and is determined by the length of tubing required. Failure to follow these recommendations could lead to pressure between generator and application.*

- |                   |   |
|-------------------|---|
| < 10 meters:      | Use 6/4 (6mm O/D, 4mm I/D) P.T.F.E. tubing.   |
| > 10 - 40 meters: | Use 10/8 (10mm O/D, 8mm I/D). Tubing and fittings not supplied in the fittings kit.   |
| > 40 metres:      | Please contact Peak Scientific with the relevant distance and we will calculate the flow resistance and the tubing size required. |

A combination of 6/4 and 10/8 tubing may be used to ensure that there is no large diameter tubing within the lab (i.e. for the first 20 meters from the Generator use 10/8 and the final 10 meters to the Mass Spectrometer use 6/4 tubing). Keep the connections and bends to a minimum.

The imperial equivalents are: 6/4 = 1/4" O/D, 3/16" I/D.  
10/8 = 3/8" O/D, 5/16" I/D.

## IMPORTANT DOCUMENTS



### Warranty Entitlement

To register your generator for your warranty entitlement, send the completed form to Peak Scientific by:

- *Email* [warranty@peakscientific.com](mailto:warranty@peakscientific.com)
- *Online* [http://www.peakscientific.com/service-and-support/warranty\\_registration](http://www.peakscientific.com/service-and-support/warranty_registration)
- *Phone* +44 (0)141 530 4185
- *Fax* +44 (0)141 812 8200

PRODUCT WARRANTY REGISTRATION	
COMPANY:	CONTACT NAME:
ADDRESS:	
	EMAIL ADDRESS:
CITY/TOWN:	<b>GENERATOR SERIAL NUMBER:</b>
POSTCODE:	
COUNTRY:	MODEL TYPE:
TELEPHONE:	<b>INSTALLATION DATE (DD/MM/YYYY):</b>

### *Important Please Note:*

You have 1 month to register your Peak Scientific product from the date of shipment.

If you wish to defer installation of your generator you must notify Peak Scientific within 1 month of the shipment date. This can be done by emailing [warranty@peakscientific.com](mailto:warranty@peakscientific.com) Once registered the warranty will be honoured for a period of 12 months after the installation date.

*For any generators that remain unregistered the warranty will begin from date of shipment.*

Thank you on behalf of Peak Scientific.

## *Normal Operation*

The Genius 3040 Gas Generator is designed specifically to minimize operator involvement. Given that the system is installed as described in earlier sections and is serviced in accordance with the specified maintenance recommendations (see [Service Requirements](#)), then it should simply be a matter of turning the Generator on when it is required.

The Generator will automatically produce the factory set flow and pressure as detailed in the [Technical Specifications](#).

## *On Demand Gas*

The system produces gas on demand. If the Mass Spectrometer is operating and requires a gas flow, the system will supply this to suit the requirements of the Mass Spectrometer. If the Mass Spectrometer requirement for gas stops, the system will also stop, once it has reached its upper set limit in the internal storage tanks. If the demand from the Mass Spectrometer starts again, the system will detect the demand for gas and will automatically start again to suit the demand.

## *Generator Cycling*

The Generator is designed for the internal compressors to cycle whilst the system is in operation and the Mass Spectrometer is demanding gas. This cycling reduces the duty (run time) on the compressors. The rate at which they cycle will be dependent on the gas required to satisfy the demand of the Mass Spectrometer. If the Mass Spectrometer demands the maximum gas flow of the system, the compressor duty will be higher, (the rest period in the compressor cycle will be shorter). If the Mass Spectrometer demand is lower than the maximum gas flow, then the duty on the compressors will decrease, (the rest period in the compressor cycle will be longer).

## *Unusual Operation*

If at any time the system begins to emit excessive noise or vibration, then it should be switched off and you should contact Peak Scientific or the Peak Partner from which the Generator has been purchased.

## Service Requirements

### Service Schedule

Service Interval	Component	Part No.	Qty.
12 months	First stage filter element	00-0031	4
	Normal acting filter	00-0032	4
	Inlet filter element	02-1054	1
	RAC filter element	00-4425	3
5000 hours	Compressor assembly	08-8316	4
Alternative to compressor assembly **	Compressor re-fit kit	06-8320	4

Table 2: Service schedule

\*\* Compressors can be re-fitted as an alternative to replacement up to a maximum of 3 times. This is a more cost effective solution, although a degree of technical expertise is required and can be time consuming. Please contact your service provider for more information.

As an alternative to purchasing the 12 month service items individually an Annual Service Kit is available as one part number. This contains all the filters required for this Generator and offers a costs saving over buying the components separately.

Purchase Interval	Component	Part No.	Qty.
12 months	Genius 3040 Annual Service Kit	08-4786	1

Table 3: Annual Service Kit

## Service Indication

The Genius 3040 counts the accumulative run-time (in hours) for the internal compressors. The Generator has the following Service Indication Stages for each system:-

### Stage 1

Once the system reaches a total of 5000 run hours, the service LED indicator (yellow) on the front of the Generator will illuminate.

This is to make the user aware that a service of the Generator is due and should be planned at the earliest convenience. The Generator will continue to operate as normal with the LED on.



*It is highly recommended that the compressors are serviced or replaced after 5000 run hours to prevent any unplanned downtime. Please contact your service provider for more information.*

## Stage 2

If the service is not completed the Generator will continue to run. Once the system reaches a total of 5250 run hours, the service LED indicator (yellow) will start to flash.

This is to make the user aware that the service of the Generator is now overdue and must be completed immediately to ensure the continuous trouble free operation of the Generator.



*Once 5250 hours is reached it is highly likely that the compressors will fail in the near future. Please contact your service provider as soon as possible to arrange a service or replacement of your compressors.*

## Service Indication Reset

Once the service has been completed the Service Indication LED can be reset in the main control PLC. This will be performed by the Peak Service Engineer or trained service representative that completes the service operation.

## Service Plans

Peak Scientific offer two service plans. The Complete Service Plan, specifically designed for Generators operated in critical environments, also includes full breakdown cover, guaranteed response times and Generator upgrades if available. Our Standard Service Plan, covering the basic needs of our Generators, features special deals on spare parts and breakdowns.

If you want to know more about our Service Plan options and how we ensure that your instrument can run uninterrupted with maximum uptime and performance, please contact us at [maintenance@peakscientific.com](mailto:maintenance@peakscientific.com)

## Cleaning

Clean the outside of the Generator only using warm soapy water and a clean damp cloth. Ensure the cloth is thoroughly rung out to remove excess fluid prior to use.



*Cleaning should only be undertaken with the power switched off and the power cord removed from the rear of the Generator.*



*Under no circumstances should any solvents or abrasive cleaning solutions be used as these can contain fumes that could be harmful to the Generator.*



## *High Duty Indication*

The Genius 3040 has a 'HIGH DUTY' indicator on the front panel. This indicator monitors the running condition of the internal compressors and illuminates when they have been operating continually for a period of 8 hours.

There are a number of extreme conditions that can affect the duty cycling of the compressors inside the Genius 3040. These extremes include very high ambient temperatures, low supply voltages, a very high flow requirement and locations of high altitude. Any one of these or a combination of a number of these extremes can be evident in a customer site.

While the Genius 3040 Generator is designed to operate fully in these conditions it should be noted that the effect of these extremes can force the compressors to run continually. The compressors are fully capable of running continually with no detrimental effect. However it should be noted that if this is the case it will increase their duty and they will reach the 5000 hour recommended service interval quicker.

## *High Duty Indication Reset*

If the compressors have been operating continually for a period of 8 hours the indicator light will be illuminated. Once the compressors return to a cycling mode the indicator light will automatically switch off. There is no manual intervention required.



## *Indication of Fault*

In most installations the 'HIGH DUTY' indicator will never be illuminated. If your Genius 3040 Generator has been operating for a significant period of time, then suddenly develops a 'HIGH DUTY' indication with no apparent change in environmental conditions or flow requirements, it may be indicative of a problem with the system such as an external leak or imminent compressor failure.

If this is the case you should contact Peak Scientific or your service provider.

*It should be noted that the Generator will continue to operate so long as it can maintain pressure during this indication.*

## Trouble Shooting

Problem	Possible Solution
The Generator will not switch on and the power switch does not illuminate.	<ul style="list-style-type: none"> <li>• Ensure power cord is plugged into the Generator and that the power socket is turned on.</li> <li>• Check the fuse in the power cord plug.</li> <li>• Contact your service provider.</li> </ul> 
The Generator will not switch on but the power switch is illuminated.	<ul style="list-style-type: none"> <li>• Disconnect power cord from the rear of the Generator. Open the left hand panel and check that both circuit breakers are turned on (switch in the up position). Reconnect power cord.</li> <li>• Contact your service provider.</li> </ul> 
Compressors are running but pressure is not building.	<ul style="list-style-type: none"> <li>• Contact your service provider.</li> </ul>
The mass spec is reporting low pressure.	<ul style="list-style-type: none"> <li>• Check pressure gauges are showing <a href="#">normal pressure</a>.</li> <li>• Contact your service provider.</li> </ul>
Yellow LED on front panel is on constantly.	<ul style="list-style-type: none"> <li>• A compressor(s) due for service. Contact your service provider.</li> <li>• Refer to <a href="#">Service Indication</a> section of this manual for further information.</li> </ul>
Yellow LED on front panel is flashing.	<ul style="list-style-type: none"> <li>• A compressor(s) is overdue for service. Contact your service provider urgently.</li> <li>• Refer to <a href="#">Service Indication</a> section of this manual for further information.</li> </ul>
Red LED on front panel is on constantly.	<ul style="list-style-type: none"> <li>• A compressor(s) has not cycled for 8 hours. Refer to the <a href="#">Run Time Alarm</a> section of this manual.</li> <li>• Ensure ambient temperatures are within the <a href="#">specification</a>.</li> <li>• Turn the Generator off and on to reset the alarm and contact your service provider.</li> </ul>
Red LED on front panel is flashing.	<ul style="list-style-type: none"> <li>• A compressor(s) has not cycled for 12 hours. Refer to the <a href="#">Run Time Alarm</a> section of this manual.</li> <li>• Ensure ambient temperatures are within the <a href="#">specification</a>.</li> <li>• Turn the Generator off and on to reset the alarm and contact your service provider.</li> </ul>
Red LED on front panel is flashing constantly and the buzzer will not stop even after power cycling the unit.	<ul style="list-style-type: none"> <li>• A compressor(s) has not cycled for 72 hours. Refer to the <a href="#">Run Time Alarm</a> section of this manual.</li> <li>• Ensure ambient temperatures are within the <a href="#">specification</a>.</li> <li>• The alarm must be reset by a qualified service technician. Contact your service provider.</li> </ul>

*Peak Scientific UK*  
Fountain Crescent  
Inchinnan Business Park  
Inchinnan  
PA4 9RE  
Scotland, UK

*Tel:* +44 (0)141 812 8100  
*Fax:* +44 (0)141 812 8200

*Peak Scientific Germany*  
Herriotstrasse 1  
60528 Frankfurt

Germany

*Tel:* +49 (0)69 677 33 490  
*Fax:* +49 (0)69 677 33 200

*Peak Scientific Taiwan*  
4F.-6, No.736  
Zhongzheng Rd.  
Zhonghe Dist.  
New Taipei City

*Tel:* +886 2 8226 2383  
*Fax:* +886 2 8226 9499

*Peak Scientific North America*  
19 Sterling Road  
Suite #1  
Billerica, MA 01862

USA

*Tel:* +1 866 647 1649  
*Fax:* +1 978 608 9503

*Peak Scientific China*  
Room 606 Building 1  
Lane 2277 Zuchongzhi  
Road  
Pudong New Area  
Shanghai  
201203, China

*Tel:* +86 21 5079 1190  
*Fax:* +86 21 5079 1191

*Peak Scientific Brazil*  
Peak Scientific Brasil  
Av. Quieroz Filho  
1700 - torre Sunny-Conj.  
602  
São Paulo SP Brasil

*Tel:* +1 866 647 1649

*Peak Scientific India*  
202, Amsri Shamira  
Old Lancer Line  
Opp. St. Mary's Degree  
College  
S.D. Road Secunderabad  
500 025, India

*Tel:* +91 40 2780 0663  
*Fax:* +91 40 2780 0663

*Peak Scientific Mexico*  
Solon 352  
Col. Los Morales Polanco  
11530 Mexico, D.F.

Mexico

*Tel:* +1 866 647 1649  
*Fax:* +1 978 608 9503

*Peak Scientific Japan K.K.*  
2-7-56, 2F Fuji Building 28  
Kita Aoyama, Minato-Ku  
Tokyo, Japan  
107-0061

*Tel:* +81 3-6864-0468

*Peak Scientific Africa*  
PO Box 478  
Somerset Mall  
7137

South Africa

*Tel:* +27 (0)2185 16542  
*Fax:* +27 (0)8654 64473

*Peak Scientific Australia*  
PO Box 65  
Belmont  
Victoria 3216

Australia

*Tel:* +61 1300 965 352

*Peak Scientific Singapore*  
3 Science Park Drive  
#03-14 The Franklin  
Science Park Drive 1  
Singapore, 118223

*Tel:* +65-6777 1966  
*Fax:* +65-6777 1978